

Quick Start Guide



Unpacking your instrument

The PFXi series is supplied with:

- PFXi series instrument (PFXi-195, PFXi-880, PFXi-950 or PFXi995)
- External power supply
- Set of 3 power leads (UK, European and United States)
- USB cable
- Ethernet cable
- Quick Start Guide
- CD ROM (containing windows® control software and manuals)
- Accessory Box

Switching on Your Instrument

Remove all packaging from your PFXi instrument; remove the silica gel pack from the sample chamber. Connect the external power supply to the 24 Volt input connector and then switch on the instrument. When the instrument is switched on, the following screen will appear.



Please make a note of the Reg. Code as this is required for the Registration of your instrument. Full details of how to register your instrument are given on the CD ROM.

CD ROM

The CD ROM supplied with your instrument contains all the User Manuals, Software and the instructions on how to perform the RCMSi calibration.

RCMSi Starter Pack

The RCMSi Starter Pack should be used within the first month of receiving the instrument. You will have to register your instrument before this calibration can be performed. Instructions on how to register your instrument and perform the RCMSi calibration are contained on the CD ROM.

The Tintometer Ltd : Lovibond House : Sun Rise Way : Amesbury : SP4 7GR : UK
T. +44 (01980) 664800: F. +44 (01980) 625412 : E. sales@tintometer.com